

## How to Use PathoCide Sanitizer/Disinfectant

### Using PathoCide on General Hard Surfaces:

This works best for phones handsets, mice, keyboards, door handles, touch plates, railings, etc. Do not spray touch screens, smart phone screens, or monitor screens.

1. Spray PathoCide onto a paper towel until wet
2. Wipe down hard surfaces
3. Let air dry

### Using Pathocide on Horizontal Hard Surfaces:

This works best for tables, counters, desks, etc.

1. Spray PathoCide onto surface
2. Wipe down with a paper towel over entire surface
3. Let air dry

### Key Factors:

- DO NOT spray chemicals directly on skin. Spray away from your face.
- Gloves are not required to use PathoCide. However, if you notice sensitization, gloves are recommended.
- Wash your hands thoroughly for 20 seconds with soap and water after using PathoCide.